


如何安全 使用排风柜 (2)

NO. 007



1. 不要在故障排风柜内进行操作
2. 在安装和拆卸设备除外，其他任何时刻不要将头部伸入排风柜内
3. 不得将排风柜台面作为化学试剂或废弃物存放场所，不要堵塞导流板槽缝
4. 排风柜内禁止使用明火
5. 减少（避免）在排风柜视窗前的人流和快速走动，不要放置风扇或带有风扇的设备
6. 避免在排风柜内使用电插线板
7. 不要将通风柜安装在门窗或空调出回风口附近

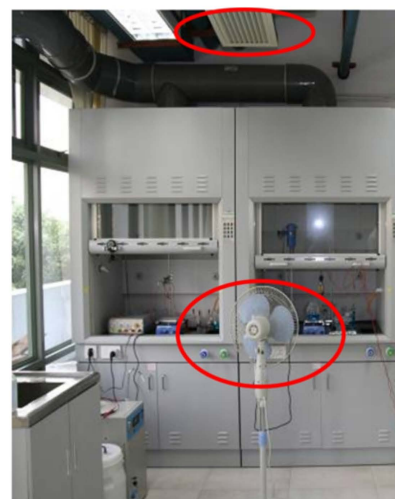

如果排风柜中有泄漏、溢出或起火...

1. 关上视窗（在安全情况）
2. 对于化学品溢出（没有火灾）事故，启动通风柜紧急通风模式（如果通风柜是有这配置）
3. 有火灾时，关上视窗，不要启动紧急通风模式
4. 联系实验室负责人和 HSE 部门求助

How to Safely Use a Fume Hood (2)



NO. 007

1. Avoid operations in faulty FM (fume hood)
2. Don't put your head in the FM unless installing or removing equipment
3. Don't store chemicals or waste on the work surface; Don't block the grooves of the baffles
4. No open flames in the FM
5. Reduce (avoid) movement near the FM, and avoid placing fan or fan-equipped devices
6. No extension cords in the FM
7. Don't place FM near doors, windows, or air vents



If a FM has a leak, spill, or fire...

1. Close the sash (safely)
2. For spills without fire, use emergency ventilation (if has)
3. For fires, close sash, and don't use ventilation
4. Contact lab supervisor and HSE for help