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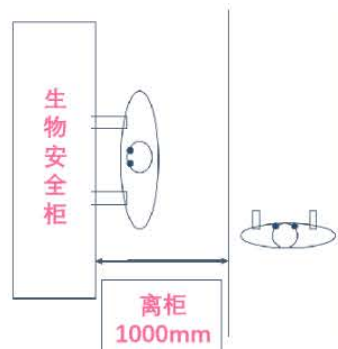
NO. 016

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

# Guidelines for Use of Biosafety Cabinets (BSC)


NO. 016

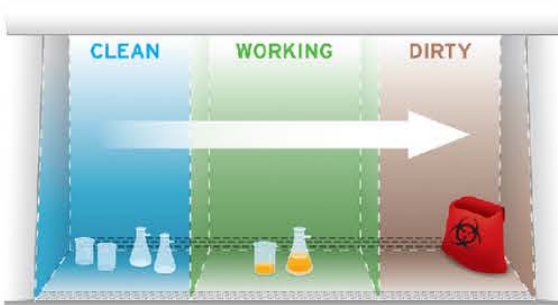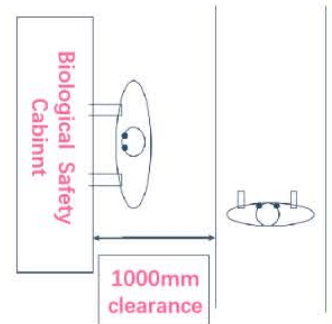- Allow the blower to run for 10~15 min before work, 5~10 min before turning off.
- Minimize disturbances to airflow curtain. Avoid rapid movements, minimize foot traffic around hood (Any pedestrian traffic routes should be at least 1000 mm from the front of the BSC). Move in a controlled and steady manner.
- Always set the sash to correct opening height (160mm~250mm for Class I, II BSC) when working in the cabinet.
- UV lamps are not substitute for good cleaning practice. Decontaminate work zone with sanitizer after use.
- DO NOT use for flammable or explosive materials as far as possible. The combustible vapor could enter the blower motor coil, and spark could cause explosion.
- When cleaning the BSC, use an alcohol-soaked towel or alcohol spray bottle to minimize the concentration of volatile solvents in the circulating air in the BSC.
- DO NOT work in the cabinet when the UV light is on. Before turning on the UV light, always make sure that the sash is fully closed to protect the user from UV radiation.
- Wear complete Personal Protective Equipment (gloves, lab coats, masks, medical caps, shoe covers, goggles, etc.).
- The cabinet with exhaust duct is required when using vaporizing toxic chemicals.
- Annual BSC Certification ensures cabinet airflow is within safe limits and the leakage rate is qualified.
- Work as deep into the work zone as possible to avoid blocking of the front or back grille (perform work at least 4 inches in front of front grille and use elbow rest). When the grille is blocked, a localized containment failure can happen that may affect the operator or contaminate the samples.

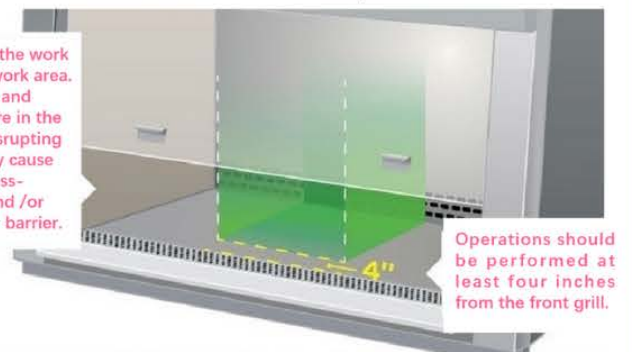The middle third of the work surface is the ideal work area. Place materials and equipment with care in the cabinet to avoid disrupting airflow, which may cause turbulence cross-contamination, and/or disruption of the air barrier.

- Dispose pipette tips and other biological waste into a medical waste garbage bag placed inside the cabinet, not outside, to prevent spreading contaminants.
- Always work "clean to dirty". Segregate sterile zone on one side, work area on middle, and waste zone on opposite side.
- DO NOT place unnecessary items inside the cabinet. Always disinfect items put into the cabinet before taking it out.
- NOT suggest to use bleach when cleaning the BSC, which can cause rust.
- DO NOT place large or multiple magnets on the BSC body. Strong magnets must not be attached to the BSC because it can disrupt the sash window sensor.
- It is strongly recommended not to use burners inside the cabinet. The hot air can disrupt the airflow causing contamination, filter damage, and even explosion.
- DO NOT plug high current equipment on the leakage protection socket inside the BSC.
- DO NOT make cabinet exhaust too close to ceiling or supply diffuser.